From glass to laboratory high-tech

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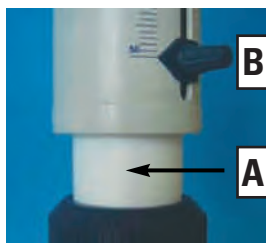
LABmax "airless"

The LABmax "airless" is the latest dispenser generation with automatic air purging.

- PTFE cylinder (A) with special ejection channel
- piston made of borosilicate glass 3.3
- adjustment screw (B) for high-precision volume adjustment

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All dispensers allow total disassembling for complete cleaning. Please follow the operating manual. All components can be replaced without problems.

Due to the constant technical improvements our products may differ from actual shown supplies.

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„Air-purging“ fully automatic

The LABmax „airless“ has an automatic air purging system. That avoids wrong values which concerns to gas bubbles in the device. Gas -lighter than liquid- is firstly leaving the device automatically through the special ejection channel. A contamination of the bottle contents is not possible because the air purging is managed through the ejection cannula.

If you do not wish air exchange between reagent and room air the LABmax is equipped with a connector which allows the simple attachment of different filters as well as calcium chloride tubes or a check valve.

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Convenient volume setting

Adjusting the volume of the LABmax is done by an adjustment screw. Just open the adjustment screw a little bit - move to the desired volume - fix the screw - thats all. Precise reproducibility of the volume in the clinical field, science and research the word is LABmax. After all - if the LABmax has to be disassembled e.g. for cleaning it does not need to be readjusted after assembling. The zero-point is part of the construction and automatically calibrated.

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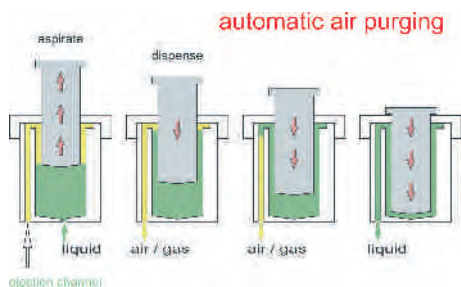
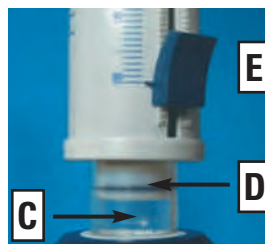
LABmax "OE"

The LABmax "OE" is the economy dispenser without automatic air-purging.

- cylinder out of borosilicate glass 3.3 (C)
- patented long life PTFE piston (D)
- fast volume adjustment with rocker switch (E) and sprocket jags (except 100 ml "OE")

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Smooth and fatigue-free dispensing

The LABmax „airless“ is made of high resistant materials and therefore high chemically and mechanically resistant and can be consequently universal (except HF). We offer the „HF“ version for hydrofluoric acid separately. The vertical valves assure a minimum of dead volume and also make effective cleaning possible. See operation manual for details.

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Safe by lock - a secure rest position

By turning the operation mode selector by 180° all liquid in the discharge tube flows back into the bottle. It is almost impossible for any leakage to occur when it is in that position.

The LABmax is completely shut down and no accidental discharge will occur by inadvertently moving the piston. This is clearly indicated by the position of the tube and the setting of the operation mode.

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LABmax „airless“

5.370.80*

LABmax airless „HF“
hydrofluoric acid
5.370.82*

LABmax "OE" without air purging system, with patented PTFE-piston, complete with accessories (see Scope of Supply)

Volume	Division steps	Tolerance	CV	Thread	Order No.
Volume	Traits de division	Tolérance	Répétabilité	Raccord à vis GL	N° de code
ml	ml	ml	ml	GL	
0.25-2.5	0.05	± 0.012	0.002	32/33	5.370.001 OE
0.5-5.0	0.10	± 0.03	0.005	32/33	5.370.002 OE
1.0-10.0	0.20	± 0.06	0.010	32/33	5.370.003 OE
2.5-25.0	0.50	± 0.15	0.025	45	5.370.004 OE
5.0-50.0	1.00	± 0.30	0.050	45	5.370.005 OE
10.0-100.0	2.00	± 0.60	0.100	45	5.370.006 OE
25.0-250.0	5.00	± 1.60	0.250	45	5.370.000 OE

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LABmax „S“ for solvents and PTFE-swelling media, with piston made of Borosilicate glass, complete with accessories (see Scope of Supply)

Volume	Division steps	Tolerance	CV	Thread	Order No.
Volume	Traits de division	Tolérance	Répétabilité	Raccord à vis GL	N° de code
ml	ml	ml	ml	GL	
1.0-10.0	0.20	± 0.06	0.010	32	5.370.013 OE
2.5-25.0	0.50	± 0.15	0.025	45	5.370.014 OE
5.0-50.0	1.00	± 0.30	0.050	45	5.370.015 OE

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LABmax „HF“ suitable for hydrofluoric acid, with patented PTFE-piston, liquids comes in contact only with PTFE and platinum iridium, complete with accessories (see Scope of Supply)

Volume	Division steps	Tolerance	CV	Thread	Order No.
Volume	Traits de division	Tolérance	Répétabilité	Raccord à vis GL	N° de code
ml	ml	ml	ml	GL	
1.0-10.0	0.20	± 0.06	0.010	32	5.370.023 OE
2.5-25.0	0.50	± 0.15	0.025	45	5.370.024 OE

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LABmax "airless" with automatic air purging system, universal use, complete with accessories (see scope of supply)

Volume	Division steps	Tolerance	CV	Thread	Order No.
Volume	Traits de division	Tolérance	Répétabilité	Raccord à vis GL	N° de code
ml	ml	ml	ml	GL	
0.25-2.5	0.05	± 0.012	0.002	32/33	5.370.801
0.5-5.0	0.10	± 0.03	0.005	32/33	5.370.802
1.0-10.0	0.20	± 0.06	0.010	32/33	5.370.803
2.5-25.0	0.50	± 0.15	0.025	45	5.370.804
5.0-50.0	1.00	± 0.30	0.050	45	5.370.805
10.0-100.0	2.00	± 0.60	0.100	45	5.370.806

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LABmax airless „HF“ suitable for hydrofluoric acid, with automatic air purging system, liquids comes in contact only with PTFE and platinum iridium, complete with accessories (see Scope of Supply)

Volume	Division steps	Tolerance	CV	Thread	Order No.
Volume	Traits de division	Tolérance	Répétabilité	Raccord à vis GL	N° de code
ml	ml	ml	ml	GL	
1.0-10.0	0.20	± 0.06	0.010	32	5.370.823
2.5-25.0	0.50	± 0.15	0.025	45	5.370.824
5.0-50.0	1.00	± 0.30	0.050	45	5.370.825

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Scope of Supply:

LABmax with operation mode selector for air-purging, complete with filling- and discharge tubes, instruction manual as well as the following thread adapters:

- 2.5/5 and 10 ml adapter A 25, A 28, A 45, S 40
- 25/50 and 100 ml adapter A 28, A 32 / 45, S 40
- conformity certification, CE
- quality certification, individual serial number
- developed and produced acc. to ISO 9001:2000

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LABmax FUTURE

- The LABmax FUTURE offers a large extent of applications for the dosage of aggressive reagents, e. g. strong acids like H_3PO_4 , H_2SO_4 , alkaline solutions like NaOH, KOH, sodium solutions and a larger number of organic solvents
- All parts coming into contact with liquid are made of following high-quality materials:
Borosilicate glass, Ceramics, Platinum-Iridium, PTFE, PFA, FEP
- Limitations of use:
Steam pressure max. 500 mbar
Viscosity max. 500 mm²/s
Temperature max. 40 °C
Density max. 2.2 g/cm³
- The LABmax FUTURE should be used in application fields of „Dispensing“, „Sampling“, „Pipetting“, „Pumping“ and is especially suitable in biochemistry use for the dispensing of culture media

LABmax FUTURE calibration inside, conformity certified, with patented PTFE-piston, complete with accessories

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Discharge tube with integrated movable tube holder

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Volume	Form	Order No.
Volume	Forme	N° de code
ml		
0.5, 2.5, 5.0, 10.0	fine pointed / fine pointue	5.370.311
0.5, 2.5, 5.0, 10.0	fine pointed / fine pointue	5.370.340
0.5, 2.5, 5.0, 10.0	standard	5.370.341

Flexible discharge tube, in spiral, made of PTFE

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Filling tube for LABmax FUTURE, made of FEP

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Filling hose for LABmax FUTURE, made of PTFE, for barrels and tanks

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Filling valve with sealing for LABmax FUTURE

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LABmax Future

5.370.114

Filling valve with sealing for LABmax FUTURE „HF“

Volume Volume ml	Order No. N° de code
10.0	5.370.660

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Discharge valve with sealing for LABmax FUTURE

Volume Volume ml	Order No. N° de code
0.5, 2.5, 5.0, 10.0	5.370.710

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Discharge valve with sealing for LABmax FUTURE „HF“

Volume Volume ml	Order No. N° de code
10.0	5.370.760

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MINISPENSOR

The bottle top dispenser for dosings in the microliter range with 2 „fix adjustable“ volumes

In medical and diagnostical laboratories a dispenser in the microliter range will be required, with which accurate volumes can be dosed with one hand only. This operation was done in the past by microliter pipettes. The top technology of witeg has fulfilled this requirement with engineering the MINISPENSOR.

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Easy to maintain

The MINISPENSOR is constructed easy to maintain. For sophisticated cleaning the upper part can be completely dismantled.

The MINISPENSOR is completely sterilizable. **Le MINISPENSOR est entièrement autoclavable**

Conformity

With the symbol  we certify the conformity with the specifications of Annex 12 of the German Weights and Measures Regulations provided the instrument is used with accessories supplied by us and in accordance to the operating manual.

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Safe volume adjustment

The MINISPENSOR will be delivered with one volume adapter, which is subdivided in 2-fix-volumes. By turning the push button the volumes can be adjusted easily and precisely.

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When turning the push-button a second volume stop will click into place which is half of the total lifting.

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Perfect handling

Unlike the conventional dispensers is the fatigue-proof handling. Uncramped, relaxed and stressless long series can be managed. With three fingers only the new developed dispenser can be manipulated. The push-button will be pressed down for dispensing with the thumb as with automatic pipettes. With the help of a spring the selected volume will be drawn back to the cylinder automatically.

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It's your money going down the drain!

Please, calculate for once your monthly consumption of pipette tips. Maybe you can exchange an expensive pipette by a MINISPENSOR.

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Safe the environment

Whenever you do not really need to work with a pipette we recommend our tip saving MINISPENSOR.

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MINISPENSOR

Fix-Volume-Bottle-Top-dispenser x 2
for accurate delivery of microliter quantities. Just screw it onto the bottle, press the plunger - thats all!

- the bottle-top micro dispenser even for hospital laboratories (glucose test, Gama GT, Chreatinine, Urea, Cholesterol, Amylase, Lypose, Iric acid, Bilirubin a.s.o.)
- for single handed operation
- due to direct displacement no sticking or jamming of the plunger
- suitable for storage in a fridge
- substitution for microliter pipettes
- special volume adapters available on request
- just select the required volume and press the plunger - thats all
- no tips required

MINISPENSOR

2x-Fixvolume-dispenser

Complete with accessories (with thread connector GL 28, 1 volume adapter and 1 bottle, both ST 28/32, suction tube and discharge tube)

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Bottles for LABmax, MINISPENSOR, TITREX
made of soda lime brown-glass, LEVASINT®-coated, screw cap made of Polypropylene (PP), space saving square shape

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Bottle for LABmax, TITREX
made of brown glass, with center screw neck and side entry filler neck for refilling the bottle without removing the LABmax or TITREX, capacity 2000 ml

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Table with 4 columns: Suitable for type / Pour type, Outer Ø / Ø ext. (mm), Inner Ø / Ø int. (mm), and Order No. / N° de code. Rows include LABmax 2.5 - 10 and LABmax 25 - 100.

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Table with 2 columns: Order No. / N° de code. Row includes 5.375.011.

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Table with 2 columns: Thread / Filetage (mm) and Order No. / N° de code. Rows include A 22, A 25, A 28, A 30, A 32 / A 28*, A 38, A 40, A 45 / A 32**, and A 45.

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Table with 2 columns: Order No. / N° de code. Row includes 5.377.222.

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Table with 2 columns: Order No. / N° de code. Row includes 5.377.300.

Calcium chloride tube, made of borosilicate glass, with plastic coating. to protect reagents in the bottle from moisture from the outside

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Order No.
N° de code
5.377.310

ST-adapters for LABmax, MINISPENSOR, TITREX, made of borosilicate glass, 1 piece each pack

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ST-adapters for
LABmax, MINISPENSOR, TITREX

WITOPED XP and WITOPED ECO

electrical pipette filling for professionals

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WITOPED XP rechargeable Pipetting for all pipettes from 0.1 to 100 ml, complete with 4 filters and recharger, unit ready for use

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WITOPED ECO the low priced pipetting help with 8 different speed levels for suction and discharging. For glass and plastic pipettes from 0.1 ml up to 100 ml. Recharger and 2 liquid stopping one-way filters included

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TITREX 2000

Burettes digital

Using the **TITREX 2000**, you will „recognize“ the changing point faster than with common glass burettes, more sensitive and, for sure less costly than with motor burettes. Hereby the dosing speed is adjusted when turning the handwheels steplessly from fast to drop by drop. The witeg burette has the advantage that the piston only without headweight is moving during the titrating process and few turnings on the wheel are sufficient to take the preselected volume. A total volume of 50 ml should be sufficient because all burette volumes can be simulated since the volume which is being sucked is determined via the display and continuously read out.

TITREX 2000

1 instrument for all volumes

**Software and interface cord
FREE OF CHARGE**

Your special advantage!

You do not need several sizes of the **TITREX 2000** - just push the button „DOS“, watch the display and with only a few turns of the handwheel you fill the burette cylinder with volumes between 0 to 50 ml - enough for several titrations.

Calibration

The permanent calibration of the **TITREX 2000** is adjusted to water at 20 °C.

For liquids of different viscosities/densities the **TITREX 2000** can be easily recalibrated by pushing the button „O/CAL“ - just follow the detailed instructions of the manual.

You can set calibration intervalls with referring reminder (recalibration timer).

Practical and fast pre-selection of volumes needed for titrations

Just push the key „DOS“ and you can pre-select the burette-volume for your titrations. By watching the display you always can control the volume of liquid, that you suck-in to the burette cylinder. Press the button „zero“ and titration can start immediately. It could not be quicker, handier and more accurate than with a glass burette and more sensitive than with a motor driven burette. The key „MEM“ (memory) helps to avoid complicated calculations.

...finally

- automatic power-cut-off after 3 minutes
- operating temperature 10-40 °C
- transport and storage temperature max. 60 °C
- a flashing indicator in the display shows, when charging of the battery is required. Running titrations remains unaffected.
- up to 1000 data sets may be stored in the memory and are available for data evaluation on PC or printers. Together with the Excel software specific evaluations are possible.
- included „Titrexsoft“-software with interface cord
- read out of operating hours

Safe and secure reposition

By turning the operation mode selector by 180° all liquid in the discharge tube flows back into the bottle. It is almost impossible for any leakage to occur when it is in that position.

The **TITREX 2000** is completely shut down and no accidental discharge will occur by inadvertently moving the piston. This is clearly indicated by the position of the tube and the setting of the operation mode.

The first Witeg burette, developed in conformity with ISO 9001 and of course conformity certified

„Calibration inside“

EN ISO 9000 ff, GLP/GMP

We have meanwhile developed a new generation of the WITEG-measuring instruments (e. g. TITREX 2000). All electronic measuring devices can justify the accuracy without any tools. Just call up the installed „CAL“ programme and the justification can be made without any problems. This device shows you continuously the new value being actualized. It's very easy and simple to work with.

All dosing and titration processes up to 1000 data can be stored non-stop and evaluated at your PC or be printed with almost any printer. Statistical data evaluation can be set up if an Excel programme is used in addition.

Titrex inclusive of Software on CD-ROM, interface cord.

TITREX 2000 digital burette, complete with filling and discharging tube, instruction manual as well as the following thread adapters: A25, A28, A32/45, S45; conformity certification. CE, single tested, quality certification, individual serial number

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5.385.200

The new witeg hand help dispenser for dosings in series in the range of 1 µl to 5000 µl, 1 WITIP filling permits up to 48 dispensing steps of 29 different volumes in intervals of 1 second without refilling and thus saves about 90 % of working time, which would have to be spent when using the traditional pipetting techniques. 1 adapter for WITIPS 25 ml and 50 ml is included to each WITOPETTE.

Technical data:
Weight: approx. 100 g
Reproducibility and accuracy:
Bidistilled water at 20-25 ± 0.5 °C constant (acc. to DIN 12 650)
WITOPETTE hand held dispenser for dosings in series from 1 µl - 5000 µl, incl. adapter for WITIPS 25 and 50 ml, for the use with WITIPS, Combitips® and Brandtips

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Table with 3 columns: Capacity, Standard pack, Order No. Rows include capacities from 0.05 to 50.0 ml and corresponding order numbers.

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Table with 2 columns: Order No., N° de code. Value: 5.385.234

WITIPS, sterile, single packed suitable for WITOPETTE and Multipette® 4780

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Table with 3 columns: Capacity, Standard pack, Order No. Rows include capacities from 0.5 to 50.0 ml and corresponding order numbers.

WITIPS, made of glass and PTFE, suitable for WITOPETTE and other dispensers - for corrosive liquids and aggressive acid p. ex triflour which are attacking the standard Polypropylene of the conventional tips

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Table with 4 columns: Capacity, Screw-thread, Standard pack, Order No. Rows include capacities 25 and 50 ml and corresponding order numbers.

WITOPET - Single channel, Fix

Microliter-pipettes, WITOPET FIX
with tip ejector, suitable for all brands of disposable tips, optimum shaft length, ergonomic design for nonfatquing operation even over longer periods, conformity certified

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WITOPET - variable Microliterpipette

Microliter-pipettes, WITOPET variable,
suitable for all brands of pipette tips, ergonomic design to fit the hand, comfortable operation overcome the fatigue and strain of repetitive pipetting, conformity certified

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
Accessories / Accessoires

Pipette support WITOPET. Suitable for al sizes of WITOPET for 7 pipettes

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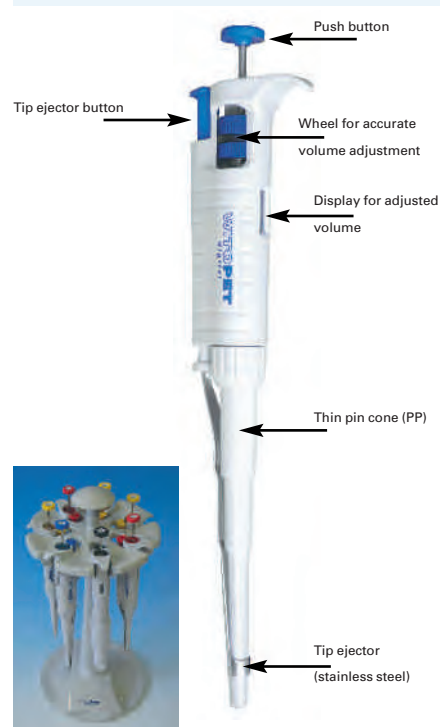WITOPET - variable economy

Microliter-pipettes, WITOPET variable economy

- Lightweight ergonomic design
- Digital display clearly reads volume setting
- The range of pipettor covers a volume range from 0.1 µl to 5000 µl
- easy to calibrate and maintenance

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WITOPET EP Digital

Microliter-pipettes, WITOPET Digital, with battery and charger console
conformity certified

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8-channel electronic pipettes, WITOPET Digital, with battery and charger console
conformity certified

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8 and 12 channel pipettes

WITOPET-Multi, variable adjustable multichannel pipette designed for comfort over long term repetitive pipetting. 8 or 12 identical volumes for titration plates.

3 volume ranges - 0.5 - 10 µl, 5 - 50 µl and 50 - 300 µl -. Highest possible accuracy and precision. Each pipette is individually tested and delivered with a test certificate.

- Designed for operator's comfort over long term repetitive pipetting
- Easy readable digital volume display, volume ranges colour coded on the push button
- Volumes steplessly adjustable
- The tip ejector can be blocked to avoid mis-application when using the instrument over large series

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5.403.1**

5.403.2**

Microliter pipettes, WITOPET-MINI, the most competitive way of quick pipetting

- with quality certificate per lot (individual certificate on request)
- Minimumsize / maximum accuracy with overall length of 130 mm only, gives the mini-pipette the complete work-feeling of a pipette
- Two stage plunger ensures last drop dispensing
- Smart tip cone design allows perfect fit for 200 µl as well as 20 µl tips
- Fully autoclavable
- Colour code
- 6 fix volumina from 5 µl to 100 µl

Volume	A	CV	Colour code	Order No.
Volume	Précision	Répétabilité	Code-couleur	N° de code
µl	≤ ± %	≤ ± %		
5	4.0	2.5	●	5.405.005
10	3.2	1.4	●	5.405.010
20	3.0	0.5	●	5.405.020
25	1.5	1.1	●	5.405.025
50	1.5	0.4	●	5.405.050
100	1.5	0.3	●	5.405.100

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WITOPETTOR, the positive displacement microliter pipette
Where air-displacement microliter pipettes have their limits, the WITOPETTOR has its field of application:

- pipetting of high viscous media, such as oils resins and fats up to viscosity of 50.000 mm²/s
- for pipetting of medias of high density up to 13.6 g/cm³ such as mercury, sulphuric acid etc.
- for pipetting of medias with foam-forming tendency
- for pipetting of liquids with high vapour pressure

WITOPETTOR variable, conformity certified, the positive lock prevents accidental volume changes

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WITOPETTOR fix, conformity certified

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Calibrated capillaries

for use with WITOPETTOR Fix and Variable, with colour code. Fire polished. Accuracy of +/-0.5%. Pack of 100 pieces

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5.406.205

Microliter-pipette tips

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Microliter-pipette Tips 0.1 to 200µl made of high quality polypropylene (PP)

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
Crystal-G Tips (0.1-10µl), short pointed tips available in bags, in racks sterile or not sterile or in Multibox Refill.


For use with Micropipettes from: Gilson Pipetman and Labmate

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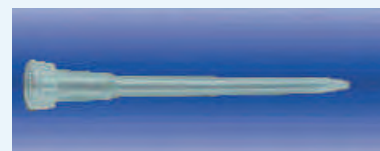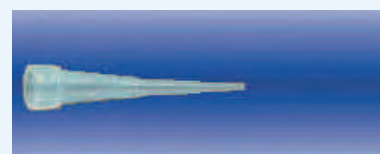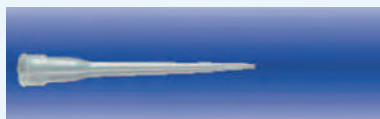
5.407.01*

5.407.02*

XL Tips colourless (0.1-10µl), long pointed tips available in bags and in racks sterile or not sterile. For use with Micropipettes from: Gilson Pipetman and Labmate

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
Yellow Tips (1-200 µl), universal long tips available in bags, in racks sterile or not sterile or in Multibox Refill.

For use with Micropipettes from: Witeg – Biohit – Brand Transferpette – Eppendorf Research or Reference – Finn/Labsystems – Gilson Pipetman – Labmate and Socorex

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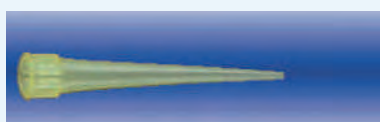Yellow Tips - G (1-200 µl), available in bags, in racks sterile or not sterile or in Multibox Refill. For use with Micropipettes from: Gilson Pipetman and Labmate

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Colourless Tips (1-200µl), universal tips available in bags, in racks sterile or not sterile or in Multibox Refill. For use with Micropipettes from: Witeg – Biohit – Brand Transferpette – Eppendorf Research or Reference – Finn/Labsystems – Gilson Pipetman – Labmate and Socorex

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Colourless Tips (1-200µl), available in bags, in racks sterile or not sterile or in Multibox Refill. For use with Micropipettes from: Labmate and Socorex

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Microliter-pipette Tips 1 to 1000µl made of high quality polypropylene (PP)

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Multi-Channel Tips (1-250µl), universal graduated tips available in bags, in racks sterile or not sterile or in Multibox Refill. For use with Multi-Channel Micropipettes from: Witeg – Biohit – Brand Transferpette – Eppendorf Research or Reference – Finn/Labsystems – Gilson Pipetman and Labmate

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Multi-Channel Tips (1-300µl), colourless universal graduated tips available in bags, in racks sterile or not sterile or in Multibox Refill. For use with Multi-Channel Micropipettes from: Witeg – Biohit – Brand Transferpette – Eppendorf Research or Reference – Finn/Labsystems – Gilson Pipetman and Socorex

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Blue Tips (200-1000 µl), universal tips available in bags, in racks sterile or not sterile or in Multibox Refill. For use with Micropipettes from: Witeg – Biohit – Brand Transferpette – Eppendorf Research or Reference – Finn/Labsystems – Gilson Pipetman – Labmate and Socorex

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Colourless Tips (200-1000 µl), universal tips available in bags, in racks sterile or not sterile or in Multibox Refill. For use with Micropipettes from: Witeg – Biohit – Brand Transferpette – Eppendorf Research or Reference – Finn/Labsystems – Gilson Pipetman – Labmate and Socorex

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Microliter-pipette Tips 200 to 5000µl made of high quality polypropylene (PP)

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Colourless Tips (500-2500 µl), tips available in bags. For use with Micropipettes from: Eppendorf Research or Reference and Socorex

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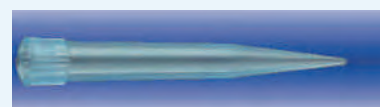
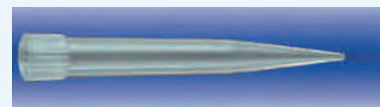
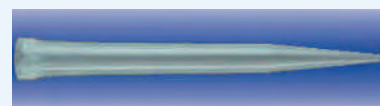
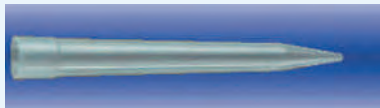
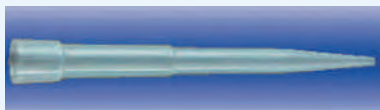
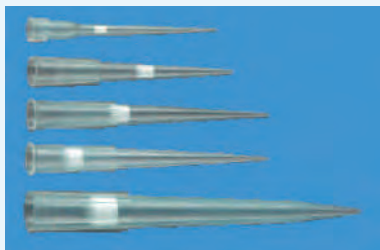
5.407.20*

5.416.000

Colourless Tips (1000-5000 µl), tips available in bags or in racks. For use with Micropipettes from: Gilson Pipetman – Labmate – Socorex

Packing Emballage	Quantity per pack Coditionnement	Order No. N° de code
Bag / <i>Sachet</i>	1 x 250	5.407.121
Rack / <i>Portoir</i>	10 x 50	5.407.122

Colourless Tips (1000-5000 µl), tips available in bags or in racks. For use with Micropipettes from: Witeg – Biohit – Brand Transferette

Packing Emballage	Quantity per pack Coditionnement	Order No. N° de code
Bag / <i>Sachet</i>	1 x 300	5.407.131
Rack / <i>Portoir</i>	10 x 50	5.407.132

Multibox plus, made of polypropylene (PP) suitable for all the multitrays, with function lid

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Multibox plus, made of polycarbonat (PC) suitable for all the multitrays, with function lid

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Microliter-pipette Filter Tips Aeroject 10 to 1000µl, made of ultra-hydrophobic PE-HD

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Microliter-pipette Filter Tips Aeroject (10-1000µl), universal, certified RNase and DNase-free and sterile packed in plastic racks with hinge lid. For use with Micropipettes from: Witeg – Biohit – Brand Transferette – Eppendorf Research or Reference – Finn/Labsystems – Gilson Pipetman – Labmate and Socorex

Volume µl	Packing Emballage	Quantity per pack Coditionnement	Order No. N° de code
10	Rack 10 x 96	960	5.407.201
20	Rack 10 x 96	960	5.407.202
100	Rack 10 x 96	960	5.407.203
200	Rack 10 x 96	960	5.407.204
1000	Rack 8 x 60	480	5.407.205

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RNase surface decontaminating agent, in spray

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Tray for reaction vessels, made of polypropylene (PP), grey, numbered from 1 - 20, for 20 reaction vessels

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